

Attach **HiScale** 50 columns to an **ÄKTA** design instrument

Instructions

Scope


Instruction on how to assemble the HiScale™ 50 column holder (28964499), and to attach the holder to an ÄKTA™ design instrument

Note: *The instruction below describes how to attach the HiScale 50 column holder to an ÄKTA avant instrument. However, the column holder can also be used on other ÄKTA design instruments.*


Instructions

Step	Action
1	<p>To assemble the HiScale 50 column holder:</p> <ul style="list-style-type: none">• Attach the column support to the column bracket with the two screws.• Insert the lobe knob through the column bracket and connect it with the screw post.

Step	Action
------	--------

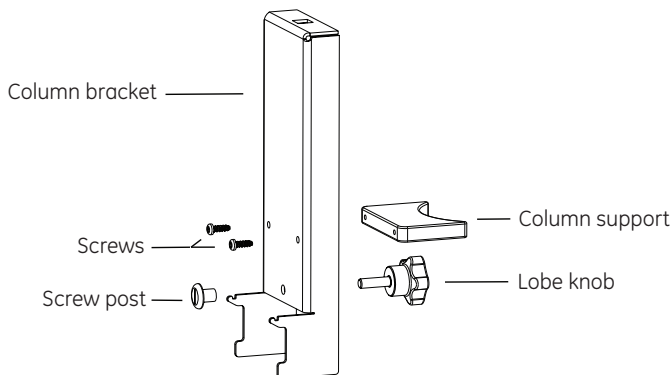


- | | |
|---|---|
| 2 | <ul style="list-style-type: none"> • Attach the column holder to a rail of the instrument by snapping-in the hooks (1) of the column bracket assembly into the rail. • Secure the column holder by turning the lobe knob (2) clock-wise. • Mount a HiScale 50 column onto the HiScale 50 column holder by placing the hook of the column in the hole on the upper surface of the holder. |
|---|---|

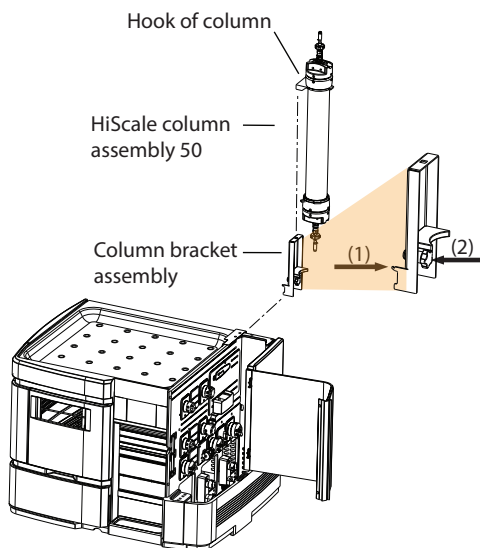

Result:

ÄKTA avant with a HiScale 50 column.

Step	Action
------	--------

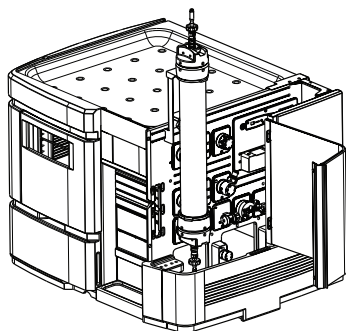



cytiva.com/akta

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

ÅKTA and HiScale are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

All other third-party trademarks are the property of their respective owners.

© 2020–2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

28976196 AB V:4 05/2021